

ProfilePrint

Accessories Box



About Us

ProfilePrint is an AI-driven ingredient quality platform powered by a patented digital food identity-as-a-service (IDaaS) technology. ProfilePrint synthesises complex molecular data of ingredient samples into digital fingerprints, rapidly predicting its identity and quality, helping agribusinesses to buy and sell better.

Backed by the world's largest food ingredient conglomerates and international venture funds, ProfilePrint is a Singapore-based company with a global footprint. For more information, please visit <https://profileprint.ai/>

If you have any question about the accessories, please contact us at support@profileprint.ai

What's in the Box?

- Specimen Dish**
x10
- Precision Specimen Spoon**
x1
- Original Specimen Spoon**
x1
- Distributor Cover**
x1
- Microfibre Cloth**
x1
- Soft Bristle Cleaning Brush**
x1

Accessories Information

This page contains information of the essential accessories for the ProfilePrint Analyser.



Standard Specimen Dish

The Standard Specimen Dish is used for solids, suspensions & colloidal specimens. Its magnetless feature is designed to seamlessly slide the dish on and off the Detachable Specimen tray. Some common ingredient examples are beans, grains and powders.

Precision Specimen Spoon

The Precision Specimen Spoon is designed to tackle more finicky ingredient types. The bowl is perfectly shaped to fit the Specimen Dish to minimise spillage during the specimen transfer from the spoon to the dish. It is suitable for the precision measurement of powders.



Original Specimen Spoon

The Original Specimen Spoon is designed to optimise speed and consistency when scooping and pouring specimens. It is suitable for high-speed scanning of solid ingredients.

Accessories Information

This page contains information of the essential accessories for the ProfilePrint Analyser.



Distributor Cover

The Distributor Cover is used to distribute and level ingredients in the Specimen Dish with a dual turn-and-compact action. The flat side of the Distributor Cover also acts as a cover for the Specimen Dish.

Microfibre Cloth

The Microfibre Cloth is used for the general cleaning of the accessories. It is recommended to be used for a final touch-up when the accessories are already relatively clean.



Soft Bristle Cleaning Brush

The Soft Bristle Cleaning Brush is used for the general cleaning of the accessories. It is recommended to be used for removing dry and solid ingredients.